

References : COMPAG100, COMPAG



100 Agarose tablets (1 tablet for 25ml gel)
Ready-to-dilute agarose tablets, instant dissolution.
Each tablet is sufficient to run two 1% gels (50 ml) or one 2% gel (25 ml)

- Add dilution buffer according to your experiments (TBE or TAE)
- Stir to dissolve the tablet
- Heat in the microwave
- After cooling (60°C) add your DNA dye (GelGreen type).
- Pour your gel

No special storage: Stable for 2 years at room temperature.

Réf : COMPAG100

20 agarose tablets (1 tablet for 25ml gel)
20 ready-to-dilute agarose tablets, instant dissolution.
Each tablet is sufficient to run two 1% gels (50 ml) or one 2% gel (25 ml)

- Add dilution buffer according to your experiments (TBE or TAE)
- Stir to dissolve the tablet
- Heat in the microwave
- After cooling (60°C) add your DNA dye (GelGreen type).
- Pour your gel

No special storage: Stable for 2 years at room temperature.

Réf : COMPAG